

Is NanoShuttle[™] biocompatible? YES!

We get asked this question all the time, and the answer is always yes. NanoShuttleTM is a nanoparticle assembly (~50 nm) consisting of gold, iron oxide, and poly-L-lysine (PLL)¹ that attaches to the plasma membrane electrostatically (50 pg/cell).

NanoShuttle™:

- Consists of biocompatible components: iron oxide and PLL are recognized as safe by the FDA^{2,3} and gold nanoparticles are in clinical trials for therapeutic use, with no indications for system toxicity⁴
- Does not bind any specific receptors, works with all cell types
- Will release off the cell over 7-8 days into the surrounding extracellular matrix, as shown by transmission electron microscopy (TEM)
- Requires magnetic forces (30 pN) only strong enough to aggregate but not harm cells
- Will not affect proliferation^{5,6}, viability⁶, metabolism^{5,7}, inflammatory⁵ or oxidative stress⁸, phenotype^{5,7,8}, and other macro cell functions
- does not cause any chromosomal abnormalities in cells, as shown by comparative genomic hybridization (CGH)

Overall, NanoShuttle™ is biocompatible and facilitates rapid 3D culture formation.

REFERENCES

- 1. Souza et al. Nat. Nanotech. (2010)
- 2. GRN No. 135, FDA. (2004)
- CFR, 21.I.A §73.200. (1994)
 ClinicalTrials.gov #NCT00848042. (2014)
- 5. Tseng et al. Tissue Eng. C. (2013)
- Daquinag et al. Tissue Eng. C. (2013)
- 7. Tseng et al. Acta Biomater. (2014)
- 8. Haisler et al. Nat. Protoc. (2013)

Spheroid Growth



Over 8 d, mCherry-tagged gioblastoma grow faster in 3D vs. $2\mathsf{D}^1$

Transmission Electron Microscopy



After 24 h (left), NanoShuttle[™] is localized with the cells, but by 8 d (right) move out of the cell and into the extracellular space¹

Proliferation, Metabolism



Neither NanoShuttle[™] nor magnetic forces have any effect on the proliferation of valvular interstitial cells (VIC, left)⁷ and 3T3 fibroblasts⁶ (right)

Viability



NanoShuttle[™] has no effect on viability, as demonstrated by live/dead staining (live = green, red = dead) on magnetically 3D bioprinted spheroids of 10,000 HepG2 hepatocellular carcinoma cells in a 384-well plate. Scale bar = 500 µm

Germany (Main office): Greiner Bio-One GmbH, info@de.gbo.com | Austria: Greiner Bio-One GmbH, office@at.gbo.com Belgium: Greiner Bio-One BVBA/SPRL, info@be.gbo.com | Brazil: Greiner Bio-One Brasil, office@br.gbo.com | China: Greiner Bio-One Suns Co. Ltd., office@cn.gbo.com

France: Greiner Bio-One SAS, infos@fr.gbo.com I Japan: Greiner Bio-One Co. Ltd., info@jp.gbo.com I Netherlands: Greiner Bio-One B.V., info@nl.gbo.com USA: Greiner Bio-One Ltd., info@uk.gbo.com WW: G

www.gbo.com/bioscience